

Resorufin Blue Cell Viability Assay Kit (Fluorometric)

LS-K241-5,000, -10,000 (5,000, 10,000 Tests) • See Storage Conditions Below



Introduction

This homogeneous assay involves simply adding a single reagent, Resorufin Blue, to the cell culture and measuring the fluorescence intensity (excitation wavelength = 530 - 570 nm, emission wavelength = 590 - 620 nm) after an incubation step. Resorufin Blue, like other resazurin-based assays such as the Alamar Blue reagent, utilizes the redox dye resazurin which is not fluorescent, but upon reduction by metabolically active cells is converted into a highly fluorescent product (resorufin). Living cells can readily reduce this non-toxic reagent and the resulting increase in fluorescence intensity can be conveniently monitored using a fluorescence spectrophotometer or plate reader. Nonviable cells have no metabolic capacity and, thus, will not reduce the dye. Therefore, the fluorescence intensity observed in this assay is a true measure of the viable cells. Resorufin Blue has been optimized for maximum sensitivity, reproducibility and long shelf-life. The homogeneous cell-based assay can be performed in multi-well plates. Resorufin Blue is compatible with all culture media and with all liquid handling systems for high-throughput screening applications in 96-well and 384-well plates. Applications include cell proliferation, cytotoxicity and apoptosis.

Key Features

- Safe. Non-radioactive assay (cf. ^3H -thymidine incorporation assay).
- Sensitive and accurate. As low as 100 cells can be accurately quantified.
- Time efficient. High-throughput assay in 96-well and 384-well plates allows simultaneous processing tens of thousands of samples per day.
- Homogeneous and convenient. A single reagent and "mix-incubate-measure" type assay. No wash and reagent transfer steps are involved.
- Robust and amenable to HTS. Z' factors of 0.6 to 0.9 are routinely observed in 96-well and 384-well plates. Can be readily automated on HTS liquid handling systems.

Applications

- Cell Proliferation: effects of cytokines, growth factor, nutrients.
- Cytotoxicity and Apoptosis: evaluation of toxic compounds, anti-cancer antibodies, toxins, environmental pollutants, etc.
- Drug Discovery: high-throughput screening for anticancer drugs.

Components

Component	K241-5000	K241-10000
	5,000 Tests	10,000 Tests
Resorufin Blue	50 mL	100 mL

Materials Not Supplied

Control Reagent: Saponin (CAS No.: 8047-15-2)

Storage

The kit is shipped at room temperature. Resorufin Blue is light sensitive and should be stored in the provided amber bottle at 4°C. The Control Reagent is stable at -20°C. Shelf life: 12 months after receipt.

FOR RESEARCH USE ONLY! Not for use in humans.

LifeSpan BioSciences, Inc. • 2401 Fourth Avenue, Suite 900, Seattle, WA 98121

www.LSBio.com • (206) 464-1554 • TechnicalSupport@LSBio.com

Resorufin Blue Cell Viability Assay Kit (Fluorometric)

LS-K241-5,000, -10,000 (5,000, 10,000 Tests) • See Storage Conditions Below



Assay Procedure

The Resorufin Blue Cell Viability assay is based on the conversion of the non-fluorescent reagent to fluorescent product by metabolically active cells. For most cells this reductive reaction takes 1 to 5 hours. The fluorescence intensity of the product is then quantified on a fluorescent microplate reader. Although most culture media contain phenol red, phenol red does not interfere with the assay. All data in Technical Notes were obtained in culture media containing phenol red.

Reagent Preparation

Important: bring reagent to room temperature before use.

Procedure Using 96-Well Plate

1. Plate and culture cells (80 μ L) in black 96-well tissue culture plates. Typical culture medium contains DMEM, 10% fetal bovine serum and antibiotics (penicillin/ strepto-mycin, gentamycin, etc), amino acids and other nutrients. Assays can be performed on either adherent cells or cells in suspension. The number of cells can vary from 100 to 80,000 per well. The volume can vary from 50 to 150 μ L, although 80 μ L is used in this protocol. In addition to the test samples, control wells of culture medium containing no cells or cells treated with a toxic reagent such as 0.1% saponin should be included.
2. Add test compounds and controls and incubate cells for the desired period of time (typically overnight). It is recommended that assays be run in duplicate or triplicate. Compounds and controls (20 μ L) can be added in phosphate buffered saline (PBS) or culture medium. The Control reagent can be conveniently reconstituted with 5 mL PBS (1% saponin).
3. Equilibrate the Resorufin Blue reagent to room temperature. Add 10 μ L (per 100 μ L of cell culture) of Resorufin Blue per well. The volume of Resorufin Blue can be adjusted depending on the volume of cell culture. Tap plate to mix cells with compounds. Incubate for 1 to 5 hours at 37°C.
4. Measure fluorescent intensity for each well on a fluorescence plate reader. If a Molecular Devices LJM Analyst is used, use the rhodamine filter sets (530nm excitation filter, 590nm emission filter and 570nm dichroic mirror).

Procedure Using 384-Well Plate

1. Plate and culture cells (40 μ L) in black 384-well tissue culture plates. The number of cells can vary from 100 to 20,000 per well. The volume can vary from 25 to 60 μ L, although 40 μ L is used in this protocol. In addition to the test samples, control wells of culture medium containing no cells or cells treated with a toxic reagent such as 0.1% saponin should be included.
2. Add test compounds and controls and incubate cells for the desired period of time. It is recommended that assays be run in duplicate or triplicate and that compounds be added in PBS or culture medium with a volume of 10 μ L.
3. Equilibrate the Resorufin Blue reagent to room temperature. Add 5 μ L Reagent (per 50 μ L of cell culture). Tap plate lightly to mix Reagent with cells. Incubate for 1 to 5 hours at 37°C.
4. Measure fluorescent intensity for each well on a fluorescence plate reader. If a Molecular Devices LJM Analyst is used, use the rhodamine filter sets (530nm excitation filter, 590nm emission filter and 570nm dichroic mirror).

FOR RESEARCH USE ONLY! Not for use in humans.

LifeSpan BioSciences, Inc. • 2401 Fourth Avenue, Suite 900, Seattle, WA 98121

www.LSBio.com • (206) 464-1554 • TechnicalSupport@LSBio.com

Resorufin Blue Cell Viability Assay Kit (Fluorometric)

LS-K241-5,000, -10,000 (5,000, 10,000 Tests) • See Storage Conditions Below



Calculations

For cell proliferation or cytotoxicity assays, the activity of a test compound can be calculated as percent change in cell number as follows,

$$\text{Activity (\%)} \text{ or Cell viability (\%)} = 100 \times (\text{F}_{\text{cmpd}} - \text{F}_0) / (\text{F}_{\text{ctrl}} - \text{F}_0)$$

where F_{cmpd} and F_{ctrl} are the average fluorescence intensities in the presence and absence (vehicle control) of the test compound and F_0 the averaged blank control fluorescence intensity.

For dose-response studies, the data are plotted against compound concentration and the EC50 for proliferation assay and IC50 for cytotoxicity assay can be determined by non-linear regression analysis using Prism or other data analysis tools.

General Considerations

Incubation time. The incubation time is dependent on the cell line. Some cell lines exhibit strong metabolic activity and, thus, require shorter incubation time than less metabolically active cell lines. The incubation time can be easily determined by reading the plate multiple times e.g. every 30 minutes after adding the Resorufin Blue reagent. In general, incubation for 1 to 5 hours is sufficient. Extensive incubation (such as >18 hours) may result in non-linear fluorescence response at high cell numbers.

Cell number. Generally the optimized Resorufin Blue reagent shows a broad range linear fluorescence response to the number of culturing cells. It is recommended to determine the number of cells per well that gives a highest signal:noise ratio. The optimal cell number can be easily determined by serial dilution of cells.

Controls. A positive control that is either cytotoxic or promotes cell proliferation can be run although it is not required. Saponin is a cytotoxic detergent. A blank control, i.e., culture medium without cells or cells containing 0.1% saponin, should be done for each assay. The blank control determines background fluorescence that must be subtracted for data analysis.

Sample Data

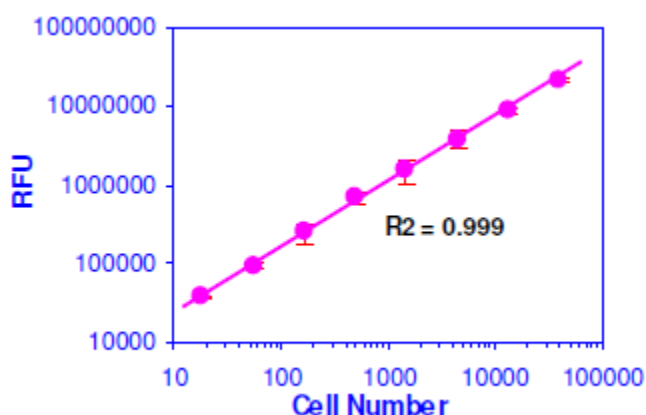


Figure 1. Linear relationship between Resorufin Blue fluorescence and cell number. HEK293 cells were serially diluted and treated with Resorufin Blue Reagent in a 384-well plate. Fluorescence intensity was linear ($R^2 = 0.999$) to the cell number. The detection limit estimated from the blank controls was 100 cells.

FOR RESEARCH USE ONLY! Not for use in humans.

LifeSpan BioSciences, Inc. • 2401 Fourth Avenue, Suite 900, Seattle, WA 98121

www.LSBio.com • (206) 464-1554 • TechnicalSupport@LSBio.com

Resorufin Blue Cell Viability Assay Kit (Fluorometric)

LS-K241-5,000, -10,000 (5,000, 10,000 Tests) • See Storage Conditions Below

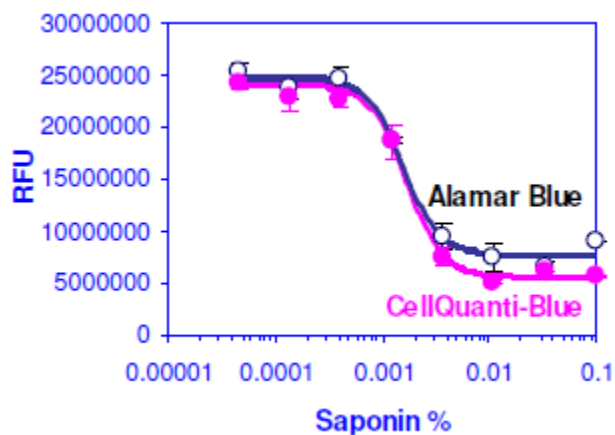


Figure 2. Dose-dependent cytotoxicity of saponin in HEK293 cells. HEK293 cells were plated at 15,000 cells per well (40 μ L). Cells were treated with varying concentrations of saponin. After culture overnight, 5 μ L Resorufin Blue reagent (solid circles) was added. Alamar Blue (open circles) was also tested for comparison. Following a three-hour incubation fluorescence intensity was measured on a Molecular Devices LJM Analyst. Data represent mean \pm SD (n = 3). The IC₅₀ values of saponin for Resorufin Blue (signal: Background = 4.8-fold) and Alamar Blue (signal: Background = 3.3-fold) were 0.0016 wt% and 0.0017 wt%, respectively.

Version: V.08.09.2018

FOR RESEARCH USE ONLY! Not for use in humans.

LifeSpan BioSciences, Inc. • 2401 Fourth Avenue, Suite 900, Seattle, WA 98121
www.LSBio.com • (206) 464-1554 • TechnicalSupport@LSBio.com