

Antioxidant Assay Kit (Colorimetric)

LS-K290-100 (100 Tests) • Store at -20°C



Introduction

An ANTIOXIDANT is a molecule capable of slowing or preventing the oxidation of other molecules. Antioxidants protect the cells from damages by reactive oxygen species which are produced in oxidation reactions in the cell. Antioxidants can be small molecules such as glutathione, vitamins, or macromolecules such as catalase, glutathione peroxidase. As oxidative stress contributes to the development of many diseases including Alzheimer's disease, Parkinson's disease, diabetes, rheumatoid arthritis and neurodegeneration, the use of antioxidants in pharmacology is intensively studied. Antioxidants are also widely used as dietary supplements and in industry as preservatives in food, cosmetics, rubber and gasoline. Simple, direct and high-throughput assays for total antioxidant capacity (TAC) find wide applications in research, food industry and drug discovery. This improved assay measures total antioxidant capacity in which Cu^{2+} is reduced by antioxidant to Cu^+ . The resulting Cu^+ specifically forms a colored complex with a dye reagent. The color intensity at 570nm is proportional to TAC in the sample.

Key Features

- Sensitive and accurate. Use 20 μL sample. Linear detection range from 1.5 to 1000 μM Trolox equivalents.
- Simple and high-throughput. The procedure involves addition of a single working reagent and incubation for 10 min. Can be readily automated as a high-throughput assay for thousands of samples per day.

Applications

- Direct Assays: serum, plasma, urine, saliva and other biological samples, food and beverages.
- Drug Discovery/Pharmacology: effects of drugs on TAC.

Components

Component	K290-100
	100 Tests
Reagent A	12 mL
Reagent B	1 mL
Standard (50 mM Trolox)	100 μL

Materials Not Supplied

Pipetting devices, centrifuge tubes, clear flat-bottom uncoated 96-well plates, plate reader capable of reading optical density at 570nm, homogenizer or sonicator etc.

Storage

The kit is shipped at room temperature. Store all components at -20 °C. Shelf life of six months after receipt.

FOR RESEARCH USE ONLY! Not for use in humans.

LifeSpan BioSciences, Inc. • 2401 Fourth Avenue, Suite 900, Seattle, WA 98121
www.LSBio.com • (206) 464-1554 • TechnicalSupport@LSBio.com

Antioxidant Assay Kit (Colorimetric)

LS-K290-100 (100 Tests) • Store at -20°C



Assay Procedure

Sample Preparation

Samples should not contain any metal chelators (e.g. EDTA) and should be clear and free of any turbidity or particles. Liquid samples (e.g. nonhemolyzed serum, plasma) can be assayed directly. Cell lysate is prepared by homogenizing or sonicating cells in ice-cold 1 x PBS and centrifugation for 10 min at 14,000 rpm to pellet any debris. Use the clear supernatant for the assay. If not assayed immediately, freeze supernatant at -80°C (stable for 1 month).

Procedure

1. Standards and Samples. Equilibrate all components to room temperature. Briefly centrifuge Reagent B and Standard before opening.

Mix 5 μL of the standard with 245 μL dH₂O (final 1 mM Trolox). Dilute standards as shown in the Table below. Transfer 20 μL standards into wells of a clear flat-bottom 96-well plate.

No	1 mM Trolox + H ₂ O	Vol (μL)	Trolox (μM)
1	100 μL + 0 μL	100	1000
2	60 μL + 40 μL	100	600
3	30 μL + 70 μL	100	300
4	0 μL + 100 μL	100	0

Transfer 20 μL of each sample into separate wells of the 96-well plate.

Note: for unknown samples, perform several dilutions to ensure that TAC is within the linear range of 1.5 to 1000 μM Trolox equivalents.

2. Assay. Prepare enough Working Reagent for Sample and Standard wells by mixing, for each assay well, 100 μL Reagent A and 8 μL Reagent B. Add 100 μL Working Reagent to all assay wells. Tap plate to mix. Incubate 10 min at room temperature.
3. Read OD_{570nm} on a plate reader.

Note: if calculated TAC is higher than 1000 μM Trolox equivalents, dilute sample in dH₂O and repeat assay. Multiply the results by the dilution factor.

Calculations

Subtract blank OD value (#4) from all standard and sample OD values. Plot the $\Delta\text{OD}_{570\text{nm}}$ against standard concentrations and determine the slope of the standard curve. Calculate the Total Antioxidant Capacity (TAC) of Sample,

$$\text{TAC} = \frac{\text{OD}_{\text{SAMPLE}} - \text{OD}_{\text{BLANK}}}{\text{Slope } (\mu\text{M}^{-1})} \times n \quad (\mu\text{M Trolox Equivalents})$$

OD_{SAMPLE} and OD_{BLANK} are the OD_{570nm} values of the sample and H₂O blank (standard #4). *n* is the sample dilution factor.

FOR RESEARCH USE ONLY! Not for use in humans.


LifeSpan BioSciences, Inc. • 2401 Fourth Avenue, Suite 900, Seattle, WA 98121

www.LSBio.com • (206) 464-1554 • TechnicalSupport@LSBio.com

Antioxidant Assay Kit (Colorimetric)

LS-K290-100 (100 Tests) • Store at -20°C

Sample Data

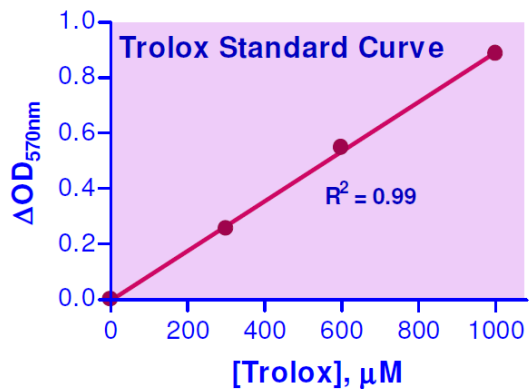

Version: V.08.09.2018

FOR RESEARCH USE ONLY! Not for use in humans.

LifeSpan BioSciences, Inc. • 2401 Fourth Avenue, Suite 900, Seattle, WA 98121

www.LSBio.com • (206) 464-1554 • TechnicalSupport@LSBio.com